


CryoSavers™

Ready-to-use Cryogenic Vials
with Brucella Broth & Cryopreservative

HARDY
DIAGNOSTICS
A Culture of Service™

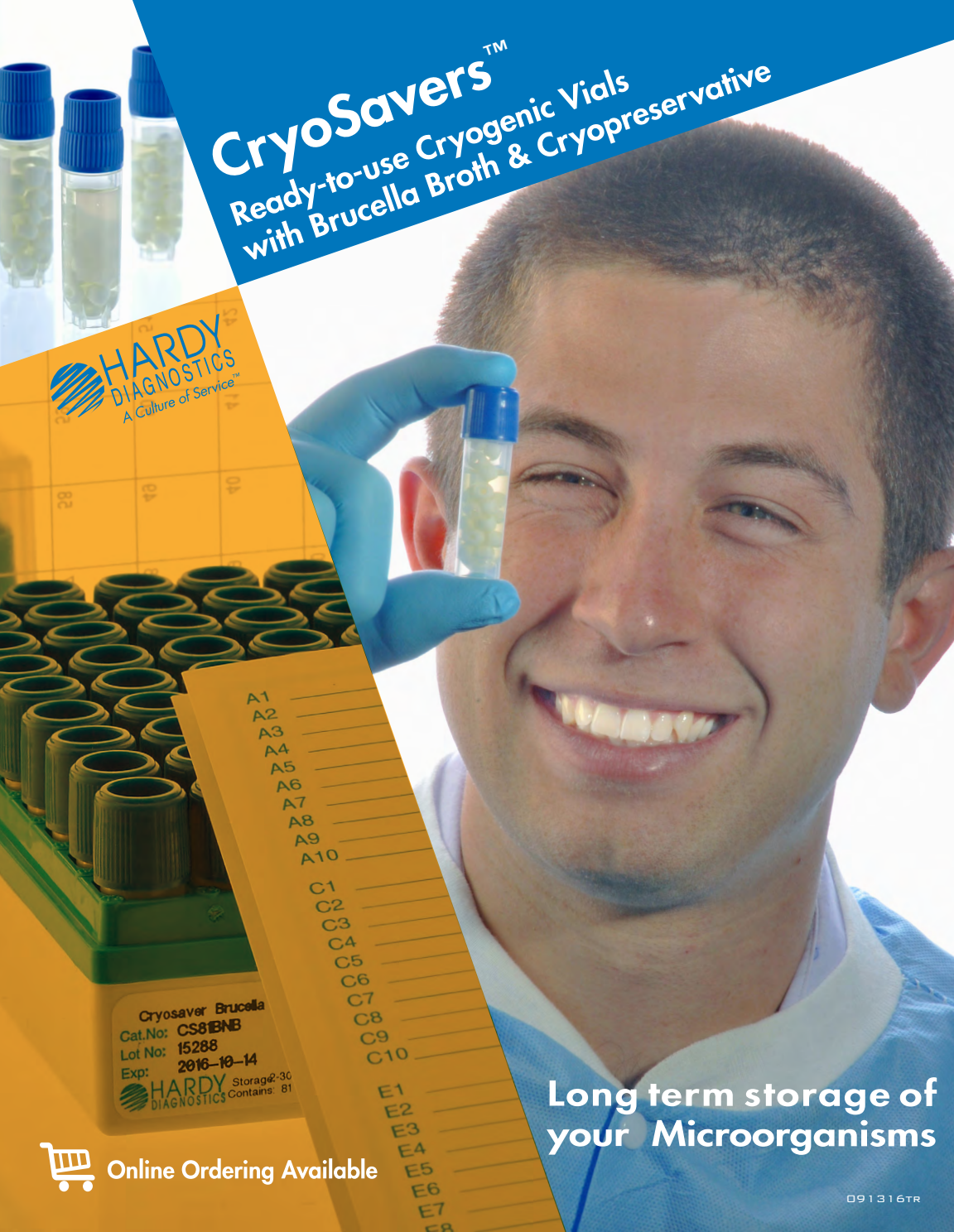

A1 _____
A2 _____
A3 _____
A4 _____
A5 _____
A6 _____
A7 _____
A8 _____
A9 _____
A10 _____
C1 _____
C2 _____
C3 _____
C4 _____
C5 _____
C6 _____
C7 _____
C8 _____
C9 _____
C10 _____
E1 _____
E2 _____
E3 _____
E4 _____
E5 _____
E6 _____
E7 _____
E8 _____

Cryosaver Brucella
Cat.No: CS81BNB
Lot No: 15288
Exp: 2016-10-14
Storage: 30
Contains: 81

Long term storage of
your Microorganisms

 Online Ordering Available



Features & Benefits

Hardy Diagnostics CryoSavers™ offers a **simple, safe,** and **convenient** way to store organisms at ultra low temperatures for future use.

Easy-to-use

Caps are ergonomically designed to feature large ridges for easy opening.

Formula Options

Available formulas include Skim Milk or Brucella Broth for the purpose of growth and long term viability of your organism.

Convenient

Ready-to-use media with cryo-preserved saves you time and expense of preparing the media yourself.

Beads

The Brucella Broth formula is available with or without plastic beads. Beads are used to roll across solid media with an inoculation needle (see "accessories").

Color Coding

Polypropylene cap inserts are also available in 11 different colors that may be placed into the top of each cap.

Accessories Available

To suit your preference, available accessories include a vial-locking workstation, spatula for scraping frozen material, and inoculation needles.

Easy Labeling

Each vial has a blank white area for written identification. Additionally, each box comes with a moisture-proof card to record the contents of individual vials.



Box contains moisture-proof log card



100/Plastic Box Cat. no. CS100BNB
81/Plastic Nalgene Box Cat. no. CS81BNB
Workstation, securely holds 40 vials, Cat. no. 802501

CryoSavers™ Vials

Type of Media	Beads	Cap Color	Cat. no.
Brucella Broth with 10% Glycerol, 1.5ml fill in a 2ml Cryovial, 100/box	No	Blue	CS100BNB
Brucella Broth with 10% Glycerol, 1.5ml fill in a 2ml Cryovial, 81/Nalgene box	No	Blue	CS81BNB
Brucella Broth with 10% Glycerol, 1.5ml fill in a 2ml Cryovial, 100/box	Yes	Blue	CS100B
Brucella Broth with Glycerol, 1.5ml fill in a 2ml Cryovial, 81/Nalgene box	Yes	Blue	CS81B
Skim Milk with Glycerol, 1.5ml fill in a 2ml Cryovial 100/box	No	Blue	CSMG100

Cap Inserts, for color coding, 500/box

Cap Insert Color	Cat. no.	Cap Insert Color	Cat. no.
White	T3121	Gray	T3129
Blue	T3122	Lilac	T31210
Red	T3123	Orange	T31211
Green	T3124	Violet	T31213
Yellow	T3125	Pink	T31214
Tan	T3128	Assorted	T3127

CryoSavers™ Product Accessories

Product Description	Cat. no.
Workstation, securely holds 40 vials, blue, 4"x 8"x 1", each	802501
Mini-inoculation needle, disposable, 1.15mm Diameter, 15cm length, blue, 25/pack, 1,000/box	COPMIN



1430 West McCoy Lane

Santa Maria, CA 93455

www.HardyDiagnostics.com | Sales@HardyDiagnostics.com

Distribution Centers: California | Washington | Utah | Arizona
Ohio | Texas | New York | Florida | North Carolina