

EnviroTrans™

ENVIRONMENTAL SWAB RINSE KIT

Swab is attached
to the cap

*A ready-to-use swab and tube pre-filled
with your choice of five solutions to
fit your specific needs!*

Collect and transport environmental samples with efficiency and ease. Simply remove the cap and attached Dacron® swab, collect the sample, recap, and send to the lab. Each EnviroTrans™ tube contains broth or buffer (see listing). The entire unit is terminally sterilized. Excellent for detection or enumeration of microorganisms from environmental surfaces or equipment.

Dacron® Tip

Available Solutions:

- **Neutralizing Buffer** for the detection of microorganisms found on surfaces. Neutralizes chlorine, iodine and quaternary ammonium compounds. Contains no growth factors, therefore is recommended for enumeration methods.
2ml, 20/box, Cat. no. SRK20
5ml, 20/box, Cat. no. SRK15
- **Lethen Broth** for the detection of microorganisms found on surfaces. Neutralizes phenols, formalin, and ethanol. May be used with 3M Petrifilm™. Contains some growth factors.
5ml, 20/box, Cat. no. SRK25
- **0.85% Saline** for the detection of microorganisms found on surfaces. Osmotically neutral transport solution. Contains no neutralizing agents or growth factors, therefore is recommended for enumeration methods.
0.5ml, 20/box, Cat. no. SRK05
2ml, 20/box, Cat. no. SRK30
5ml, 20/box, Cat. no. SRK35
- **D/E Neutralizing Broth** for the detection of microorganisms found on surfaces. Neutralizes a broad spectrum of antiseptics and disinfectants, such as quaternary ammonia, iodine, chlorine, phenolics, formalin, mercurials, glutaraldehyde, ethanol, and hexachlorophene. Detects organisms remaining after sanitation treatment. Contains some growth factors.
5ml, 20/box, Cat. no. SRK45
- **Neutralizing Saline** for the detection of microorganisms found on surfaces. Neutralizes a broad spectrum of antiseptics and disinfectants such as; quaternary ammonia, iodine, chlorine, phenolics, formalin, mercurials, glutaraldehyde, ethanol, and hexachlorophene. Detects microorganisms remaining after sanitation treatment. Contains no growth factors, therefore is recommended for enumeration methods.
5ml, 20/box, Cat. no. SRK55



*5ml D/E
Neutralizing Broth*

Self-standing Vial

Tween is a registered trademark of ICI Americas, Inc. Petrifilm is a trademark of the 3M Corporation. Dacron is a registered trademark of DuPont, Corp.

**Call for a
Free Evaluation
Sample!**

Other fills and formulas are available. Call our Technical Support department at 800.266.2222 ext. 5598 for more information.

**Call us for free evaluation samples
800.266.2222 ext. 5604**

**Sales@HardyDiagnostics.com
www.HardyDiagnostics.com**

**HARDY**
DIAGNOSTICS
A Culture of Service™



EnviroTrans™

ENVIRONMENTAL SWAB RINSE KIT



I would like more information on this product.

Please fill in your information below and fax the completed card to us at 805.346.2760.

Name _____

Title _____

Company _____

Address _____

City _____

State/Zip _____

Phone _____

Email _____

For more information on EnviroTrans™ or any other Hardy Diagnostic Product:

Call toll free at 800.266.2222 ext. 5604

Email us at Sales@HardyDiagnostics.com

Visit us online at www.HardyDiagnostics.com

A Culture of Service™



HARDY
DIAGNOSTICS

1430 West McCoy Lane
Santa Maria, CA 93455